




Overview: model class **G-ULTIMATE-90**

Safe and approved storage of flammable or toxic gases in working areas – for up to 4 x 50-litre gas cylinders



Maximum fire resistance

- G-ULTIMATE-90 cabinets offer maximum safety when storing gas cylinders – 90 minutes fire resistance

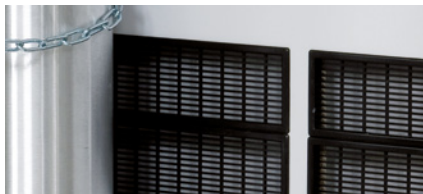

Optimal ventilation that complies with the standard

- tested and certified in accordance with the tightened GS regulations
- complete and even ventilation in the whole interior of the cabinet
- pressure drop values of all models < 150 Pa



Space-saving talent

- unique on the market: gas cylinder cabinet with a 60 cm width for the storage of two 50 litre steel gas cylinders or two 40 litre aluminium gas cylinders with a diameter of 230 mm

Further features of this model class



Locking

- doors lockable with profile cylinder lock (integrated in the door handle)
- cabinets can be integrated by the customer in an existing master key system



Pipe lead-throughs

- large area for flexible lead-through possibilities on the top of the cabinet
- no additional insulation of piping necessary



Comfortable adjusting possibilities

- integrated adjusting aids in the cabinet base ensure a fast and effective compensation for uneven floor of up to 10 mm

Possible interior equipment



- **standard interior equipment** complete with mounting rails, cylinder retainer and matching tension belts and rolling ramp



- **comfort interior equipment** complete with mounting rails, cylinder retainer and matching tension belts and rolling ramp with pneumatic damper



- **lateral cylinder retainer** for two 10-litre cylinders, alternatively in height-adjustable version (no tools required)



- **shelves** in different versions, can also be installed at an angle for the storage of small gas cylinders



Safety storage cabinets for the storage of pressurised gas cylinders in working areas according to EN 14470-2 and TRGS 510



- **fire resistance 90 minutes** (type G90), type-tested in accordance with EN 14470-2



- **GS-tested** (EN 14470-1, EN 14727, document EK5/AK4 09-10, ProdSG)
- **50,000** wear-free opening and closing actions
- test basis: all cabinets of this model group correspond to the requirements of the Equipment and Product Safety Act (§7 Para. 1) with respect to the guarantee of safety and health



- **excellent products with special manufacturer warranty**
- extended **manufacturer warranty of up to 5 years** in combination with the asecos BASICPlus tariff



G-LINE



G-ULTIMATE-90 ◀ 140 cm ▶



Gas cylinder cabinet G-ULTIMATE-90 model G90.205.140

in Light grey (RAL 7035), interior equipment with 1 x standard interior equipment for gas cylinders (sheet steel galvanised and powder coated)

Order No. 30641-001-30642

Safe and approved storage of flammable or toxic gases in working areas – for up to 4 x 50-litre gas cylinders

Function / construction:

- **no unauthorised use:** doors lockable with profile cylinder (integration in an existing locking system possible)
- **easy alignment:** adjusting aids to compensate for uneven floor
- **easy mounting of connecting pipes and gas fittings:** large interior height (1890 mm), many lead-through possibilities on the top of the cabinet
- **ventilation:** integrated air ducts ready for connection (DN 75) to a technical exhaust, even ventilation inside the cabinet
- **tested and certified:** according to the stricter GS principles


Available equipment

- comfort interior equipment – complete with mounting rails, rolling ramp (L = 350 mm) with pneumatic damper, cylinder retainer and matching tension belts
- standard interior equipment – complete with mounting rails, rolling ramp, cylinder retainer and matching tension belts
- cylinder retainer – variable installation in the cabinet, flexibly adjustable in depth, across the entire width of the cabinet
- lateral cylinder retainer – for two 10-litre cylinders, alternatively in height adjustable version (no tools required)



Lateral cylinder retainer

For two 10-litre cylinders, alternatively in height-adjustable version (no tools required)

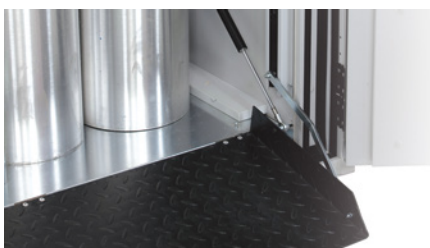

Comfort interior equipment

Extremely flat rolling angle, optimum slip resistance, folding out almost weightlessly



Complete ventilation

In the interior from the bottom to the top




Cylinder lock integrated in the door handle

The cabinet can be integrated into a locking system





Gas cylinder trolley

Direct placement of gas cylinders into asecos gas cylinder cabinets

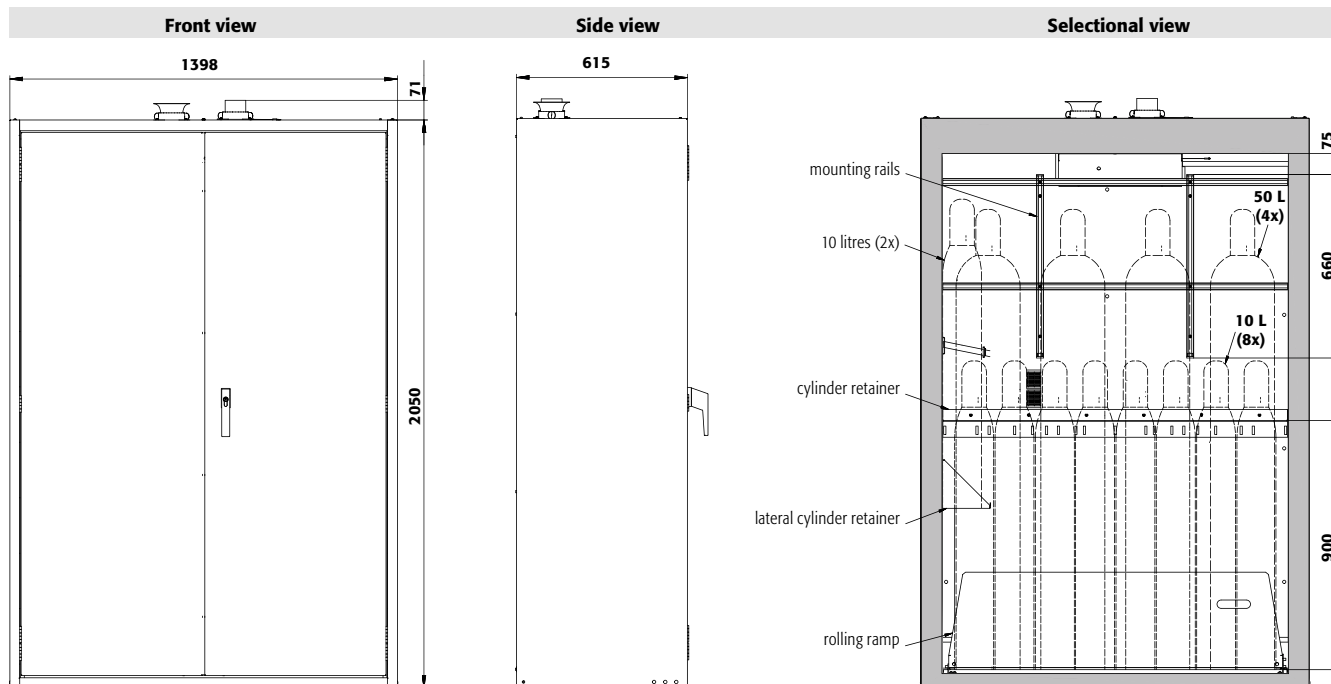**Technical data**

External dimensions W x D x H	mm	1398 x 615 x 2050
Internal dimensions W x D x H	mm	1245 x 400 x 1858
Weight without interior equip.	kg	690
Extraction air	DN	75
Air change 10 times	m ³ /h	9
Diff. pressure (air change 10 times)	Pa	5
Diff. pressure (air change 10 times EK5/AK4)	Pa	5
Air change 120 times	m ³ /h	111
Diff. pressure (air change 120 times)	Pa	54
Diff. pressure (air change 120 times EK5/AK4)	Pa	143
Maximum load	kg	600
Distributed load	kg/m ²	571
50-litre gas cylinders	piece	4
Lead-through possibilities		72

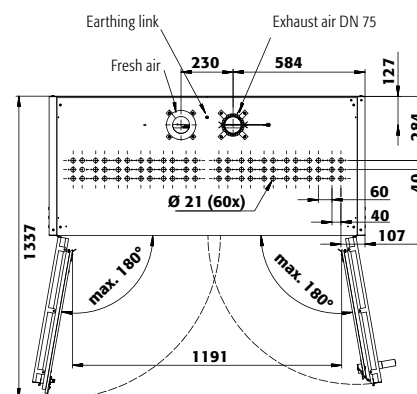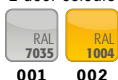G90.205.140**Top view**2 door colours are available **at no extra cost**:

001 002

Order example - how to order correctly

30641 — 001 — 30642

	Model	Information / equipment	Order No.		
	G90.205.140	without interior equipment	30641		
	Body colour	Door colour	Order No.		
	Light grey RAL 7035	Light grey RAL 7035	001 		
	Warning yellow RAL 1004	Warning yellow RAL 1004	002 		
	Interior equipment packages (can only be ordered in combination with the cabinet)	Material	Order No.		
	1x standard interior equipment for gas cylinders	sheet steel galvanised and powder coated	30642 		
	1x comfort Interior equipment for gas cylinders	sheet steel galvanised and powder coated	30643 		
	Interior equipment	Material	W x D x H (mm)	 1)  2)  3)	Order No.
	shelf	sheet steel powder-coated RAL 7035	1235 x 241 x 25*	75.00	30644 
	shelf, inclined for storage of small bottles	sheet steel powder-coated RAL 7035	1236 x 377 x 25*	75.00	30399 
	grid, as storage level	steel galvanised	1233 x 280 x 30	180.00	30645 
	cylinder retainer on the sidewall	sheet steel powder-coated RAL 7035	350 x 175 x 175		30647 
	cylinder retainer on the sidewall, height adjustable without tools	sheet steel powder-coated RAL 7035	350 x 175 x 175		30648 
	securing chain, in place of tension belts	steel galvanised			30646 
	document holder, DIN A5	polystyrene			30793 
	Accessories				Order No.
	Ermeto coupling				5981 
	tension belt, additional				13181 
	cylinder retainer, profile as a cylinder retainer incl. 2 tension belts and mounting material flexible mounting in height at the rear wall				30650 
	preparation for lead-through with female thread				30649 
	gas cylinder trolley				26294 

* W x D usable storage space / 1) load capacity (kg), 2) capacity (litres), 3) maximum container volume inside the cabinet (litres)



G-ULTIMATE-90 ◀ 120 cm ▶



Gas cylinder cabinet G-ULTIMATE-90 model G90.205.120

in Light grey (RAL 7035), interior equipment with 1 x comfort interior equipment for gas cylinders (sheet steel galvanised and powder coated)

Order No. 30657-001-30659 

Safe and approved storage of flammable or toxic gases in working areas – for up to 4 x 50-litre gas cylinders

Function / construction:

- **no unauthorised use:** doors lockable with profile cylinder (integration in an existing locking system possible)
- **easy alignment:** adjusting aids to compensate for uneven floor
- **easy mounting of connecting pipes and gas fittings:** large interior height (1890 mm), many lead-through possibilities on the top of the cabinet
- **ventilation:** integrated air ducts ready for connection (DN 75) to a technical exhaust, even ventilation inside the cabinet
- **tested and certified:** according to the stricter GS principles


Available equipment

- comfort interior equipment – complete with mounting rails, rolling ramp (L = 350 mm) with pneumatic damper, cylinder retainer and matching tension belts
- standard interior equipment – complete with mounting rails, rolling ramp, cylinder retainer and matching tension belts
- cylinder retainer – variable installation in the cabinet, flexibly adjustable in depth, across the entire width of the cabinet
- lateral cylinder retainer – for two 10-litre cylinders, alternatively in height adjustable version (no tools required)




Lateral cylinder retainer

For two 10-litre cylinders, alternatively in height-adjustable version (no tools required)

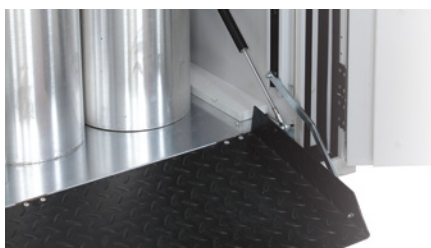

Comfort interior equipment

Extremely flat rolling angle, optimum slip resistance, folding out almost weightlessly

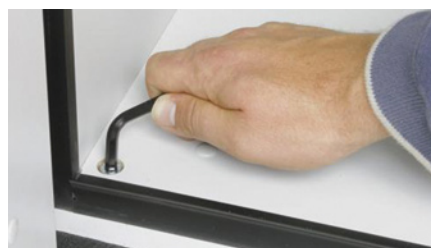

Adjusting aids

To compensate for uneven floor and an easy alignment



Complete ventilation

In the interior from the bottom to the top




Cylinder lock integrated in the door handle

The cabinet can be integrated into a locking system




Large interior height

Easy and effective mounting of the connecting pipes and gas equipment in the cabinet

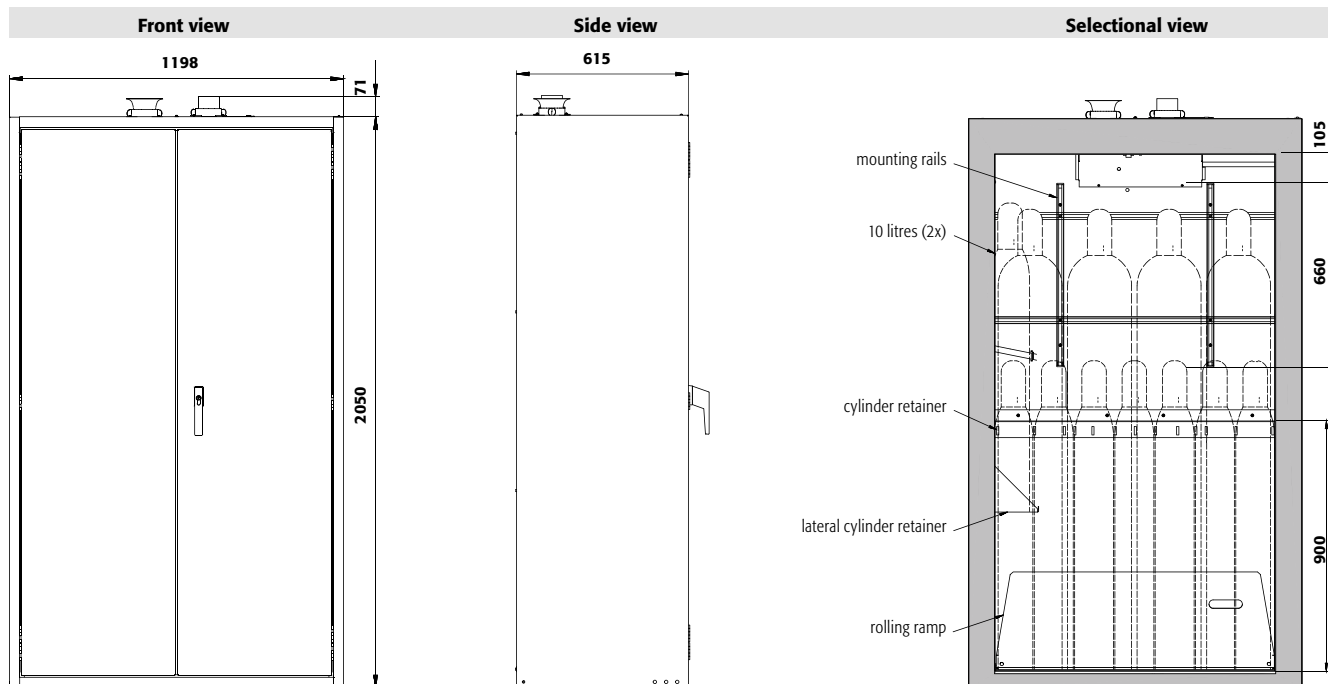**Technical data**

External dimensions W x D x H	mm	1198 x 615 x 2050
Internal dimensions W x D x H	mm	1045 x 400 x 1858
Weight without interior equip.	kg	610
Extraction air	DN	75
Air change 10 times	m ³ /h	8
Diff. pressure (air change 10 times)	Pa	5
Diff. pressure (air change 10 times EK5/AK4)	Pa	5
Air change 120 times	m ³ /h	93
Diff. pressure (air change 120 times)	Pa	49
Diff. pressure (air change 120 times EK5/AK4)	Pa	103
Maximum load	kg	600
Distributed load	kg/m ²	625
50-litre gas cylinders	piece	4
Lead-through possibilities		60

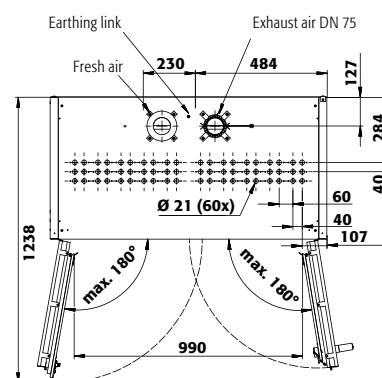G90.205.120**Top view**2 door colours are available **at no extra cost**:**Order example - how to order correctly**

	Model	Information / equipment	Order No.				
	G90.205.120	without interior equipment	30657				
	Body colour	Door colour	Order No.				
	Light grey RAL 7035	Light grey RAL 7035	001 				
	Warning yellow RAL 1004	Warning yellow RAL 1004	002 				
	Interior equipment packages (can only be ordered in combination with the cabinet)	Material	Order No.				
	1x standard interior equipment for gas cylinders	sheet steel galvanised and powder coated	30658 				
	1x comfort Interior equipment for gas cylinders	sheet steel galvanised and powder coated	30659 				
	Interior equipment	Material	W x D x H (mm)	 1)	 2)	 3)	Order No.
	shelf	sheet steel powder-coated RAL 7035	1035 x 241 x 25*	75.00			30660 
	shelf, inclined for storage of small bottles	sheet steel powder-coated RAL 7035	1040 x 432 x 30*	75.00			30771 
	grid, as storage level	steel galvanised	1033 x 280 x 30	180.00			30661 
	cylinder retainer on the sidewall	sheet steel powder-coated RAL 7035	350 x 175 x 175				30647 
	cylinder retainer on the sidewall, height adjustable without tools	sheet steel powder-coated RAL 7035	350 x 175 x 175				30648 
	securing chain	steel galvanised					30662 
	document holder, DIN A5	polystyrene					30793 
	Accessories						Order No.
	Ermeto coupling						5981 
	tension belt, additional						13181 
	cylinder retainer, profile as a cylinder retainer incl. 2 tension belts and mounting material flexible mounting in height at the rear wall						30650 
	preparation for lead-through with female thread						30649 
	gas cylinder trolley						26294 

* W x D usable storage space / 1) load capacity (kg), 2) capacity (litres), 3) maximum container volume inside the cabinet (litres)



G-ULTIMATE-90 ◀ 090 cm ▶



Gas cylinder cabinet G-ULTIMATE-90 model G90.205.090

in Light grey (RAL 7035), interior equipment with 1 x standard interior equipment for gas cylinders (sheet steel galvanised and powder coated)

Order No. 30667-001-30668 4


Safe and approved storage of flammable or toxic gases in working areas – for up to 3 x 50-litre gas cylinders

Function / construction:

- **no unauthorised use:** doors lockable with profile cylinder (integration in an existing locking system possible)
- **easy alignment:** adjusting aids to compensate for uneven floor
- **easy mounting of connecting pipes and gas fittings:** large interior height (1890 mm), many lead-through possibilities on the top of the cabinet
- **ventilation:** integrated air ducts ready for connection (DN 75) to a technical exhaust, even ventilation inside the cabinet
- **tested and certified:** according to the stricter GS principles

Available equipment

- comfort interior equipment – complete with mounting rails, rolling ramp (L = 350 mm) with pneumatic damper, cylinder retainer and matching tension belts
- standard interior equipment – complete with mounting rails, rolling ramp, cylinder retainer and matching tension belts
- cylinder retainer – variable installation in the cabinet, flexibly adjustable in depth, across the entire width of the cabinet
- lateral cylinder retainer – for two 10-litre cylinders, alternatively in height adjustable version (no tools required)

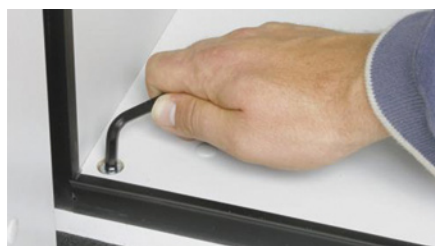

Adjusting aids

To compensate for uneven floor and an easy alignment



Lateral cylinder retainer

For two 10-litre cylinders, alternatively in height-adjustable version (no tools required)



Large interior height

Easy and effective mounting of the connecting pipes and gas equipment in the cabinet




Complete ventilation

In the interior from the bottom to the top





Gas cylinder trolley

Direct placement of gas cylinders into asecos gas cylinder cabinets

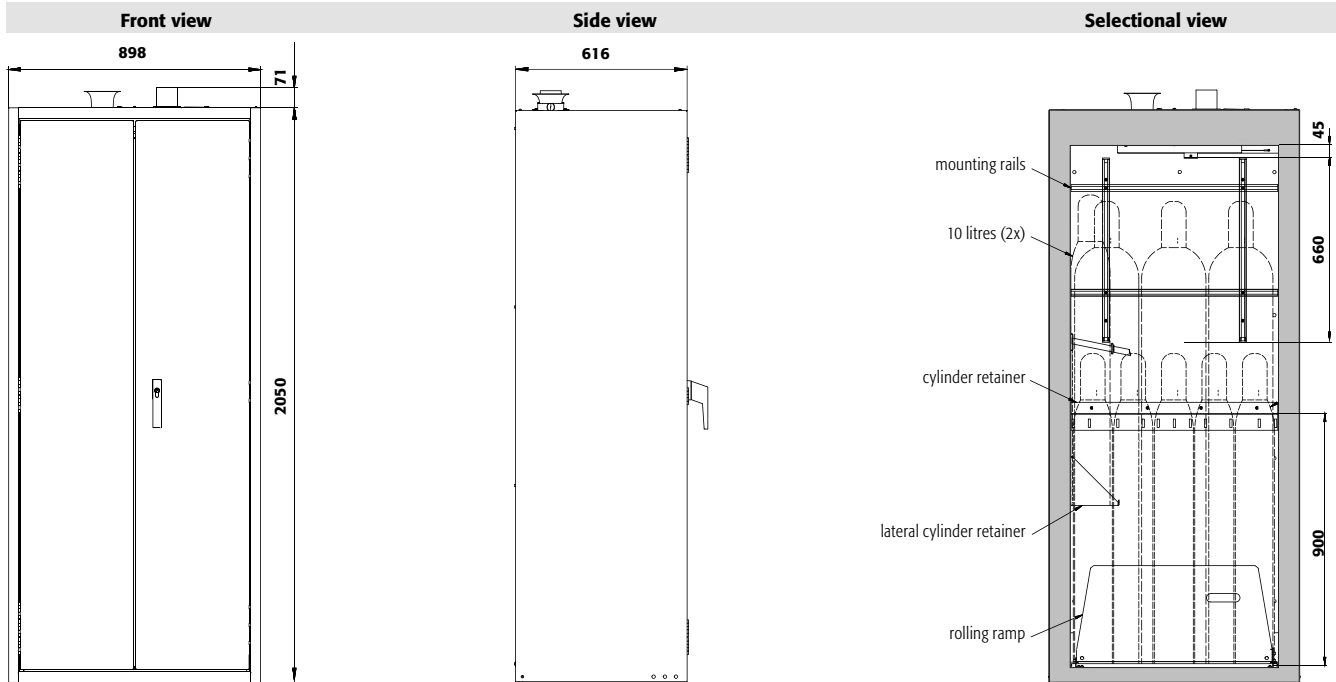**Technical data**

External dimensions W x D x H	mm	898 x 615 x 2050
Internal dimensions W x D x H	mm	745 x 425 x 1858
Weight without interior equip.	kg	490
Extraction air	DN	75
Air change 10 times	m ³ /h	6
Diff. pressure (air change 10 times)	Pa	5
Diff. pressure (air change 10 times EK5/AK4)	Pa	5
Air change 120 times	m ³ /h	66
Diff. pressure (air change 120 times)	Pa	42
Diff. pressure (air change 120 times EK5/AK4)	Pa	147
Maximum load	kg	600
Distributed load	kg/m ²	752
50-litre gas cylinders	piece	3
Lead-through possibilities		36

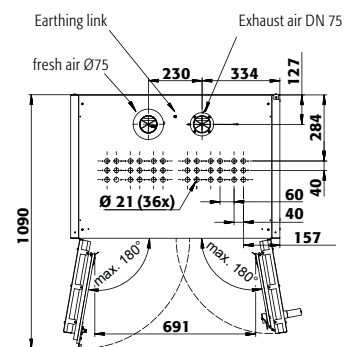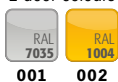G90.205.090**Top view**2 door colours are available **at no extra cost**:**Order example - how to order correctly**

	Model	Information / equipment	Order No.				
	G90.205.090	without interior equipment	30667				
	Body colour	Door colour	Order No.				
	Light grey RAL 7035	Light grey RAL 7035	001 				
	Warning yellow RAL 1004	Warning yellow RAL 1004	002 				
	Interior equipment packages (can only be ordered in combination with the cabinet)	Material	Order No.				
	1x standard interior equipment for gas cylinders	sheet steel galvanised and powder coated	30668 				
	1x comfort interior equipment for gas cylinders	sheet steel galvanised and powder coated	30669 				
	Interior equipment	Material	W x D x H (mm)	 1)	 2)	 3)	Order No.
	shelf	sheet steel powder-coated RAL 7035	735 x 241 x 25*	75.00			30670 
	shelf, inclined for storage of small bottles	sheet steel powder-coated RAL 7035	740 x 432 x 30*	75.00			30787 
	grid, as storage level	steel galvanised	733 x 280 x 30	180.00			30671 
	cylinder retainer on the sidewall	sheet steel powder-coated RAL 7035	350 x 175 x 175				30647 
	cylinder retainer on the sidewall, height adjustable without tools	sheet steel powder-coated RAL 7035	350 x 175 x 175				30648 
	securing chain, in place of tension belts	steel galvanised					30672 
	document holder, DIN A5	polystyrene					30793 
	Accessories						Order No.
	Ermeto coupling						5981 
	tension belt, additional						13181 
	cylinder retainer, profile as a cylinder retainer incl. 2 tension belts and mounting material flexible mounting in height at the rear wall						30650 
	preparation for lead-through with female thread						30649 
	gas cylinder trolley						26294 

* W x D usable storage space / 1) load capacity (kg), 2) capacity (litres), 3) maximum container volume inside the cabinet (litres)



G-ULTIMATE-90 ◀ 060 cm ▶



Safe and approved storage of flammable or toxic gases in working areas – for up to 2 x 50-litre gas cylinders

Function / construction:

- **no unauthorised use:** door lockable with profile cylinder (integration in an existing locking system possible)
- **easy alignment:** adjusting aids to compensate for uneven floor
- **easy mounting of connecting pipes and gas fittings:** large interior height (1890 mm), many lead-through possibilities on the top of the cabinet
- **ventilation:** integrated air ducts ready for connection (DN 75) to a technical exhaust, even ventilation inside the cabinet
- **tested and certified:** according to the stricter GS principles

Available equipment

- standard interior equipment – complete with mounting rails, rolling ramp, cylinder retainer and matching tension belts



Gas cylinder cabinet

G-ULTIMATE-90 model G90.205.060.2F


in Light grey (RAL 7035), interior equipment with 1 x standard interior equipment for gas cylinders (sheet steel galvanised and powder coated)

Order No. 30678-001-30682 



Lateral cylinder retainer

For two 10-litre cylinders, alternatively in height-adjustable version (no tools required)

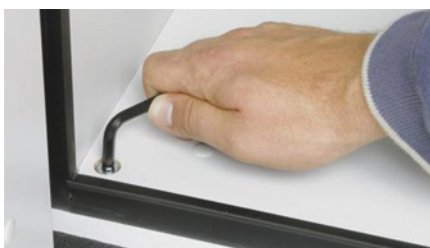

Adjusting aids

To compensate for uneven floor and an easy alignment



Cylinder lock integrated in the door handle

The cabinet can be integrated into a locking system



Large interior height

Easy and effective mounting of the connecting pipes and gas equipment in the cabinet




Gas cylinder trolley


Direct placement of gas cylinders into asecos gas cylinder cabinets




Front view

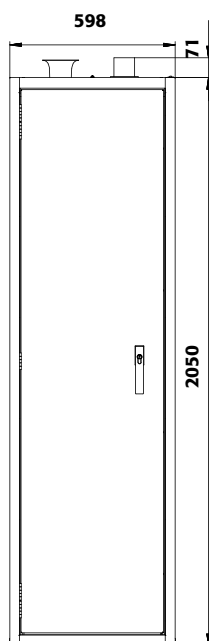

Side view

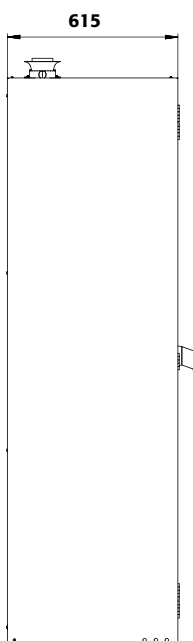

Selectional view

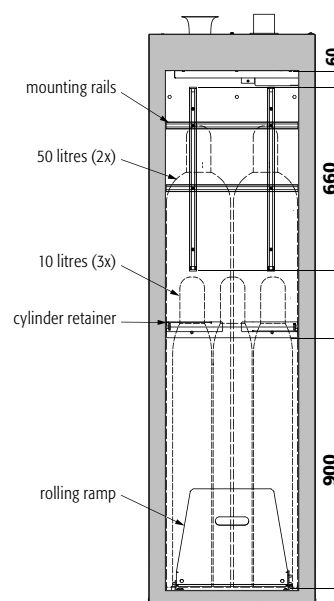


Technical data

External dimensions W x D x H	mm	598 x 615 x 2050
Internal dimensions W x D x H	mm	477 x 425 x 1858
Weight without interior equip.	kg	365
Extraction air	DN	75
Air change 10 times	m ³ /h	4
Diff. pressure (air change 10 times)	Pa	5
Diff. pressure (air change 10 times EK5/AK4)	Pa	5
Air change 120 times	m ³ /h	45
Diff. pressure (air change 120 times)	Pa	45
Diff. pressure (air change 120 times EK5/AK4)	Pa	76
Maximum load	kg	600
Distributed load	kg/m ²	999
50-litre gas cylinders	piece	2
Lead-through possibilities		24

G90.205.060.2F

Top view

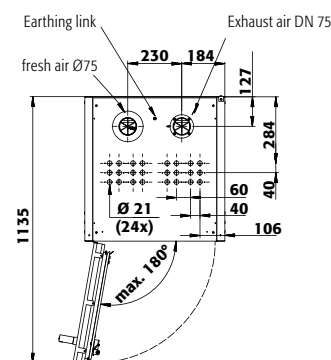

2 door colours are available **at no extra cost**:



Order example - how to order correctly



Model

G90.205.060.2F
G90.205.060.2FR

Information / equipment

without interior equipment, for 2x 50 liter cylinders, left hinged
without interior equipment, for 2x 50 liter cylinders, right hinged

Order No.

30678
30679



Body colour

Light grey RAL 7035
Warning yellow RAL 1004

Door colour

Light grey RAL 7035
Warning yellow RAL 1004

Order No.

001 **4**
002 **10**



Interior equipment packages (can only be ordered in combination with the cabinet)

1x standard interior equipment for gas cylinders

Material

sheet steel galvanised and powder coated

Order No.

30682 **4**



Accessories

Ermeto coupling
preparation for lead-through with female thread
gas cylinder trolley

Order No.

5981 **2**
30649 **2**
26294 **2**

* W x D usable storage space / 1) load capacity (kg), 2) capacity (litres), 3) maximum container volume inside the cabinet (litres)



G-ULTIMATE-90 ◀ 060 cm ▶



Gas cylinder cabinet G-ULTIMATE-90 model G90.205.060

in Light grey (RAL 7035), interior equipment with 1 x standard interior equipment for gas cylinders (sheet steel galvanised and powder coated)

Order No. 30676-001-30680

Safe and approved storage of flammable or toxic gases in working areas – for up to 1 x 50-litre gas cylinders or 2 x 10-litre gas cylinders

Function / construction:

- **no unauthorised use:** door lockable with profile cylinder (integration in an existing locking system possible)
- **easy alignment:** adjusting aids to compensate for uneven floor
- **easy mounting of connecting pipes and gas fittings:** large interior height (1890 mm), many lead-through possibilities on the top of the cabinet
- **ventilation:** integrated air ducts ready for connection (DN 75) to a technical exhaust, even ventilation inside the cabinet
- **tested and certified:** according to the stricter GS principles

Available equipment


- comfort interior equipment – complete with mounting rails, rolling ramp (L = 350 mm) with pneumatic damper, cylinder retainer and matching tension belts
- standard interior equipment – complete with mounting rails, rolling ramp, cylinder retainer and matching tension belts
- cylinder retainer – variable installation in the cabinet, flexibly adjustable in depth, across the entire width of the cabinet
- lateral cylinder retainer – for two 10-litre cylinders, alternatively in height adjustable version (no tools required)



Gas cylinder cabinet G-ULTIMATE-90 model G90.145.060

in Light grey (RAL 7035), interior equipment with 1 x standard interior equipment for gas cylinders (sheet steel galvanised and powder coated)

Order No. 30689-001-30691

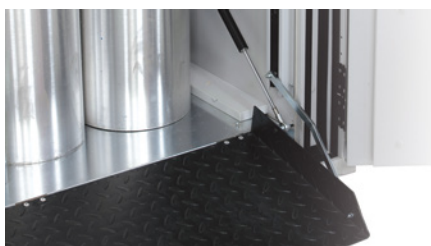

Comfort interior equipment

Extremely flat rolling angle, optimum slip resistance, folding out almost weightlessly




Cylinder lock integrated in the door handle

The cabinet can be integrated into a locking system




Gas cylinder trolley

Direct placement of gas cylinders into asecos gas cylinder cabinets

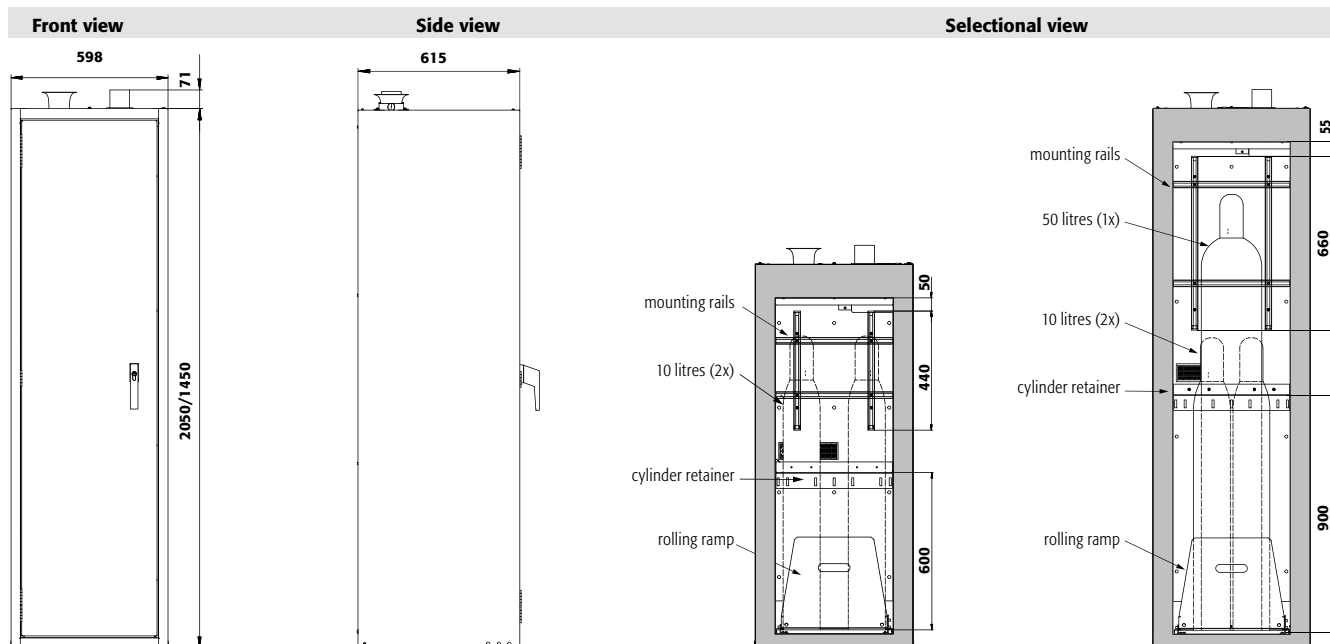**Technical data**

		G90.205.060	G90.145.060
External dimensions W x D x H	mm	598 x 615 x 2050	598 x 615 x 1450
Internal dimensions W x D x H	mm	445 x 425 x 1858	445 x 425 x 1246
Weight without interior equip.	kg	365	278
Extraction air	DN	75	75
Air change 10 times	m ³ /h	4	2
Diff. pressure (air change 10 times)	Pa	5	5
Diff. pressure (air change 10 times EK5/AK4)	Pa	5	5
Air change 120 times	m ³ /h	42	28
Diff. pressure (air change 120 times)	Pa	95	18
Diff. pressure (air change 120 times EK5/AK4)	Pa	85	42
Maximum load	kg	600	600
Distributed load	kg/m ²	999	909
50-litre gas cylinders	piece	1	
Lead-through possibilities		24	24

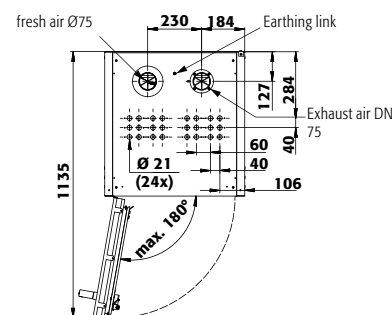Top view

Model	Information / equipment	Order No.
G90.205.060	without interior equipment, left hinged	30676
G90.145.060	without interior equipment, left hinged	30689
G90.205.060.R	without interior equipment, right hinged	30677
G90.145.060.R	without interior equipment, right hinged	30690

Body colour	Door colour	Order No.
Light grey RAL 7035	Light grey RAL 7035	001 4
Warning yellow RAL 1004	Warning yellow RAL 1004	002 10

Interior equipment packages (can only be ordered in combination with the cabinet) G90.205.060(R)	Material	Order No.
1x standard interior equipment for gas cylinders	sheet steel galvanised and powder coated	30680 4
1x comfort interior equipment for gas cylinders	sheet steel galvanised and powder coated	30681 6

Interior equipment packages (can only be ordered in combination with the cabinet) G90.145.060(R)	Material	Order No.
1x standard interior equipment for gas cylinders	sheet steel galvanised and powder coated	30691 4

Interior equipment	Material	W x D x H (mm)	1) 2) 3)	Order No.
shelf	sheet steel powder-coated RAL 7035	435 x 241 x 25*	75.00	30683 2
shelf, inclined for storage of small bottles	sheet steel powder-coated RAL 7035	440 x 432 x 30*	75.00	30788 2
cylinder retainer on the sidewall	sheet steel powder-coated RAL 7035	350 x 175 x 175		30647 2
cylinder retainer on the sidewall, height adjustable without tools	sheet steel powder-coated RAL 7035	350 x 175 x 175		30648 2


Interior equipment	Material	W x D x H (mm)	1) 2) 3)	Order No.
securing chain, in place of tension belts	steel galvanised			30684 2

Accessories	Order No.
Ermeto coupling	5981 2
tension belt, additional	13181 2
cylinder retainer, profile as a cylinder retainer incl. 2 tension belts and mounting material flexible mounting in height at the rear wall	30650 2
preparation for lead-through with female thread	30649 2
document holder, DIN A5	30793 2
gas cylinder trolley	26294 2

* W x D usable storage space / 1) load capacity (kg), 2) capacity (litres), 3) maximum container volume inside the cabinet (litres)




Model class G-ULTIMATE-90 - special solutions

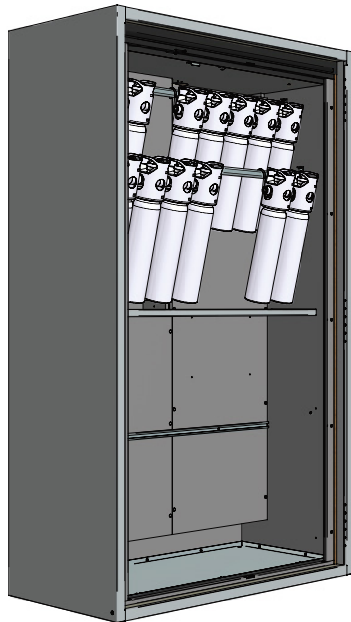


Suspended storage

of special gas cylinders, taking model G90.205.120 as an example

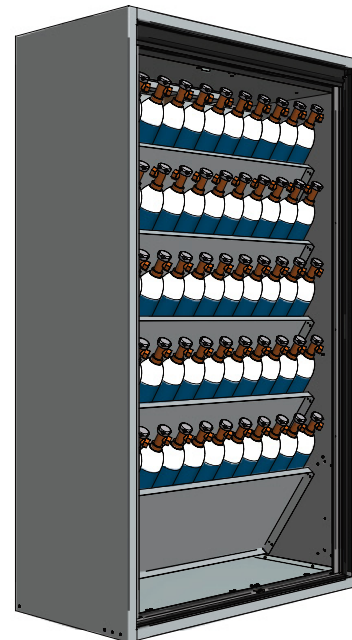


Storage of 2-litre cylinders (max. number)

taking model G90.205.120 as an example, with 5 inclined shelves
(Order No. 30771)

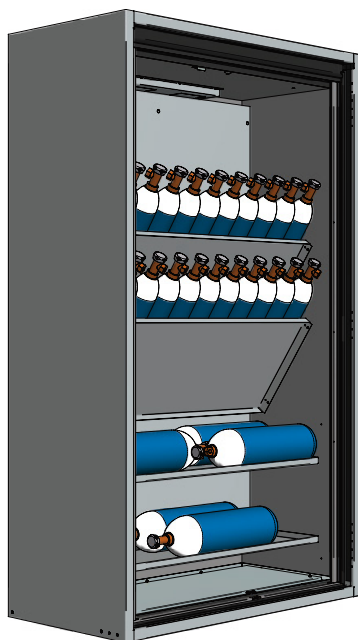

Combined storage

of 3 and 5-litre cylinders, taking model G90.205.120 as an example,
with 2 inclined shelves (Order No. 30771) and 2 shelves
(Order No. 30660)

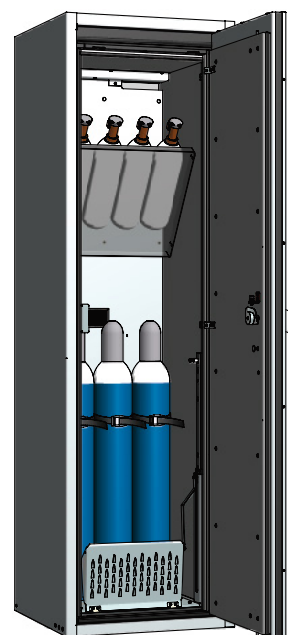

Combined storage

of 2 and 10-litre cylinders, taking model G90.205.060.R as an example,
and 1 inclined shelf with standard interior equipment (Order No. 30782)
and 1 inclined shelf (Order No. 30788)

371



WolfLabs

Pricing on any accessories shown can be found by keying the part number into the search box on our website.

The specifications listed in this brochure are subject to change by the manufacturer and therefore cannot be guaranteed to be correct. If there are aspects of the specification that must be guaranteed, please provide these to our sales team so that details can be confirmed.

www.wolflabs.co.uk

Tel : 01759 301142

Fax : 01759 301143

sales@wolflabs.co.uk

Please contact us if this literature doesn't answer all your questions.