

# Cole-Parmer®

## Cole-Parmer® Hot Plate, acid resistant, HP-200

- A completely new concept in hot plates
- PTFE construction with glass ceramic plate for exceptional resistance to chemical attack
- Ideal for acid digestions or trace metal analysis
- Separate control box connected with a PTFE coated lead



HP-200 Hot Plate

# Cole-Parmer® Hot Plate, acid resistant, HP-200



HP-200 Hot Plate



Cole-Parmer® A heated glass ceramic plate mounted in a block of pure PTFE creates a powerful hotplate which is almost impervious to chemical attack, even by concentrated acids.

When boiling acid solutions the HP-200 is unaffected by the fumes and splashes which eventually destroy conventional hotplates. The chemical inertness of the PTFE body and ceramic top plate also means that much more aggressive cleaning agents can be used. For example, if all traces of metal must be removed, this hotplate can be washed with concentrated nitric acid!

The separate temperature controller is connected to the hotplate via a 2 metre PTFE coated lead. This allows the hotplate to be located in a fume cupboard and the controller kept outside, well away from the corrosive environment. The controller is also fitted with a “Hot” warning light that will flash whenever the plate temperature of the hotplate is above 50°C and it will continue to operate when the hotplate is turned off and connected to the electricity supply.

The hotplate has a large 30x30cm square heated area so is ideal for heating either one large vessel or several smaller ones.

## Technical Specification

Specification	HP-200-AR
Plate material	Glass ceramic
Body material	PTFE
Plate dimensions, mm	300 x 300
Heated area, mm	200 x 200
Heater power, W	900
Max. plate temperature, °C	400
Hotplate dimensions, mm, (w x d x h)	320 x 360 x 60
Control unit dimensions, mm, (w x d x h)	150 x 160 x 65
Net weight, kg	11
Electrical supply	230V, 50-60Hz
IP Rating	43 (Plate) & 30 (Control)

## Ordering Information

Description	Ordering Number	Series No.	Model No.	Legacy Sku.
Cole-Parmer® Hotplate, acid resistant, including control unit	04801-08	HP-200	HP-200-AR	CP300



# WolfLabs

**Pricing on any accessories shown can be found by keying the part number into the search box on our website.**

The specifications listed in this brochure are subject to change by the manufacturer and therefore cannot be guaranteed to be correct. If there are aspects of the specification that must be guaranteed, please provide these to our sales team so that details can be confirmed.

**[www.wolflabs.co.uk](http://www.wolflabs.co.uk)**

**Tel : 01759 301142**

**Fax : 01759 301143**

**[sales@wolflabs.co.uk](mailto:sales@wolflabs.co.uk)**

Please contact us if this literature doesn't answer all your questions.